Ohaus Corporation
19A Chapin Road
Pine Brook, NJ 07058, USA
Tel: (973) 377-9000
Fax: (973) 944-7177
www.ohaus.com

Ohaus CL Series Scales
Models CL201, CL2000 and CL5000
Básculas Ohaus Serie CL
Modelos CL201, CL2000 y CL5000
Balances Ohaus série CL
Modèles CL201, CL2000 et CL5000

Operating Manual
Manual de Funcionamiento
Manuel d'utilisation

Register your scale at www.ohaus.com
under *Warranty Registration*

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© Ohaus Corporation 2004, all rights reserved.

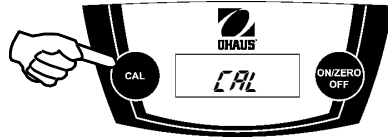
Printed in China P/N 80010609


Calibration/Calibración/Étalonnage

Refer to Optional Accessories for the correct calibration weights.

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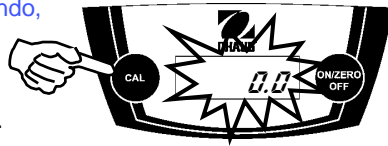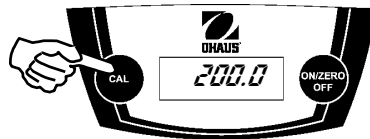
Note: If “E” is displayed a calibration error has occurred (for example, an incorrect weight was used or the pan was disturbed). Repeat the procedure.

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Install 3-AA batteries (provided), or connect the AC adapter (optional accessory) to the power input jack at the rear of the scale. Do not apply excessive force or press on the pan.

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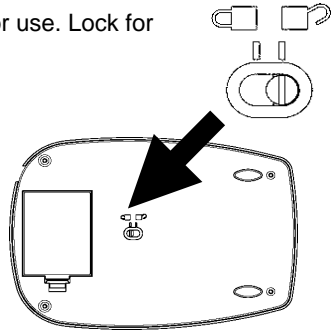
Slide the load cell lock to the unlock position for use. Lock for stacking, storing, or transport.

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
ON/ZERO/Off

Press this key to turn the scale on. This key is also used to rezero the display. Press and hold this key to turn the scale off.

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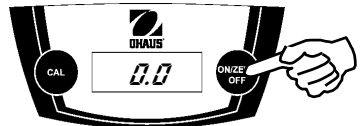Place sample on the pan and read the displayed result. The * indicator appears when the reading is stable. If using a container, press ON/ZERO/Off to tare the container before adding material.

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Specifications/Especificaciones /Spécifications

Model	CL201	CL2000	CL5000
Capacity (g)	200	2000	5000
Readability (g)	0.1	1	2
Repeatability (Std. Dev.) (g)	0.1	1	2
Linearity (g)	±0.1	±1	±2
Weighing Units	grams		
Tare range	to capacity by subtraction		
Stabilization time	3 seconds		
Optimum operating temp.	65° - 77°F/18° - 25°C		
Power requirements	3-AA alkaline batteries (included) AC adapter (optional accessory)		
Battery Life	40 hours		
Auto Shut-off	2 minutes no activity		
Overload indication	[E]		

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Ohaus products are warranted against defects in materials and workmanship for duration of the warranty. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus. This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of product registration, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages. As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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Limited Warranty/Garantía limitada /Garantie limitée